



SM1PD1B

Descriptions

The Thorlabs SM1PD1B photodiode is ideal for measuring both pulsed and CW fiber light sources, by converting the optical power to an electrical current. The detector is mounted in a convenient SM1 (Ø1.035"-40) externally threaded tube for easy mounting and integration in to existing set up. The detector also has internal RMS (Ø0.8"-36) threading for further mounting solutions. The photodiode is a *Type B (Anode Grounded)* arrangement and the pin codes for the specific package can be found in the drawing below. The photodiode anode produces a current, which is a function of the incident light power (P) and the wavelength (λ). The responsivity, $\mathcal{R}(\lambda)$, can be read from the curve Responsivity Curve (provided below) to estimate the amount of photocurrent to expect. This current can be converted to a voltage by placing a load resistor (R_L) from the photodiode anode to the circuit ground. The output voltage is derived as:

$$V_o = P \times \mathcal{R}(\lambda) \times R_L$$

The bandwidth, f_{BW} , and the rise time response, t_R , are determined from the diode capacitance, C_j , and the load resistance, R_L , as shown below. The diode capacitance can be lowered by placing a bias voltage from the photodiode cathode to the circuit ground. Ω

$$f_{BW} = \frac{1}{(2\pi)R_L C_j}, t_R = \frac{0.35}{f_{BW}}$$


Specifications



Specification		Value
Wavelength Range	λ	400 - 1100 nm
Peak Wavelength	λ_p	970 nm
Responsivity	$\mathcal{R}(\lambda)$	0.65 A/W
Active Area Diameter		94 mm ²
Rise/Fall Time ($R_L=50 \Omega$, 20 V)	t_r/t_f	40 ns
NEP, Typical (440 nm, 20 V)	W/√Hz	2.07×10^{-13}
Dark Current (5 V)	I_d	1.05 nA
Capacitance (5 V)	C_j	432 pF
Package		SM1, External Thread RMS, Internal Thread
Sensor Material		Si
Corresponding Unmounted Diode		FDS1010

Maximum Rating	
Max Bias (Reverse) Voltage	20 V
Reverse Current	10 mA
Operating Temperature	-10 to 60 °C
Storage Temperature	-20 to 70 °C

Recommended Circuit Diagram

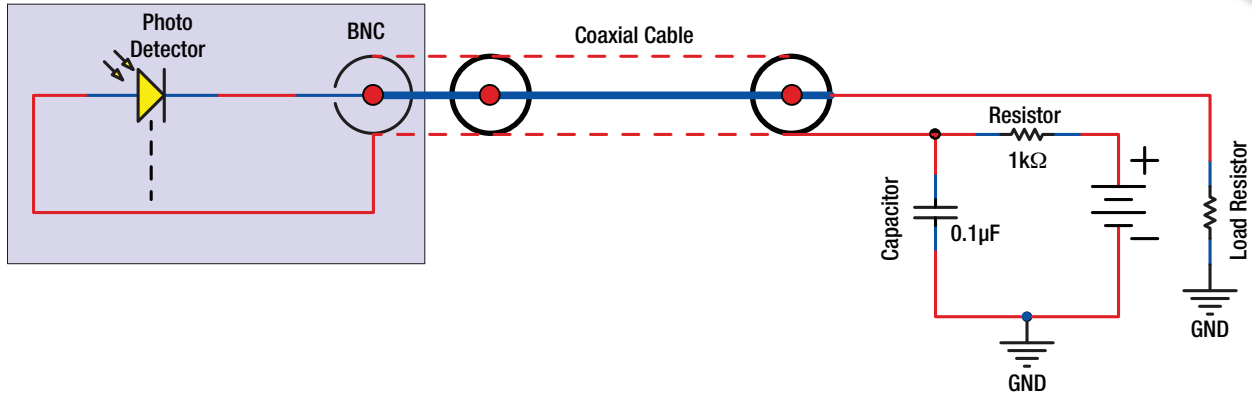


Responsivity Graph


The responsivity of a photodiode is a measure of its sensitivity to light and is defined as the ratio of the photocurrent I_p to the incident light power P at a given wavelength:

$$R_\lambda = \frac{I_p}{P}$$

In other words, it is a measure of the effectiveness of the conversion of light power into electrical current. Responsivity varies from lot to lot and with the wavelength of the incident light, applied reverse bias, and temperature. It increases slightly with applied reverse bias due to improved charge collection efficiency in the photodiode. The change in temperature increases or decreases the width of the band gap and varies inversely with the temperature.

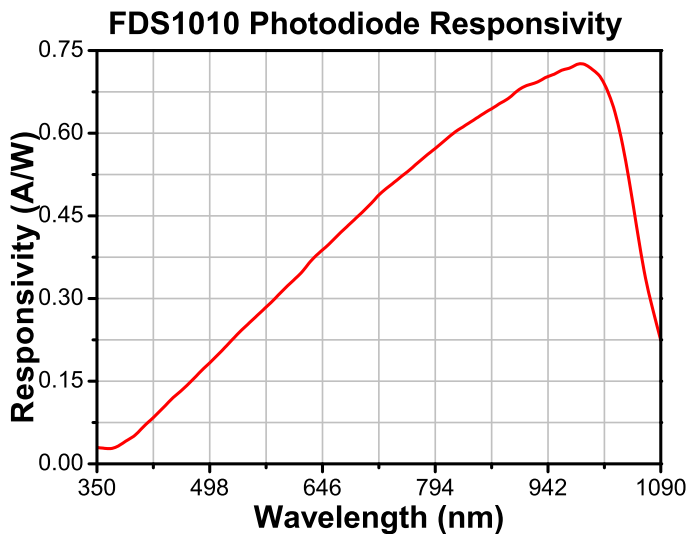

Drawing

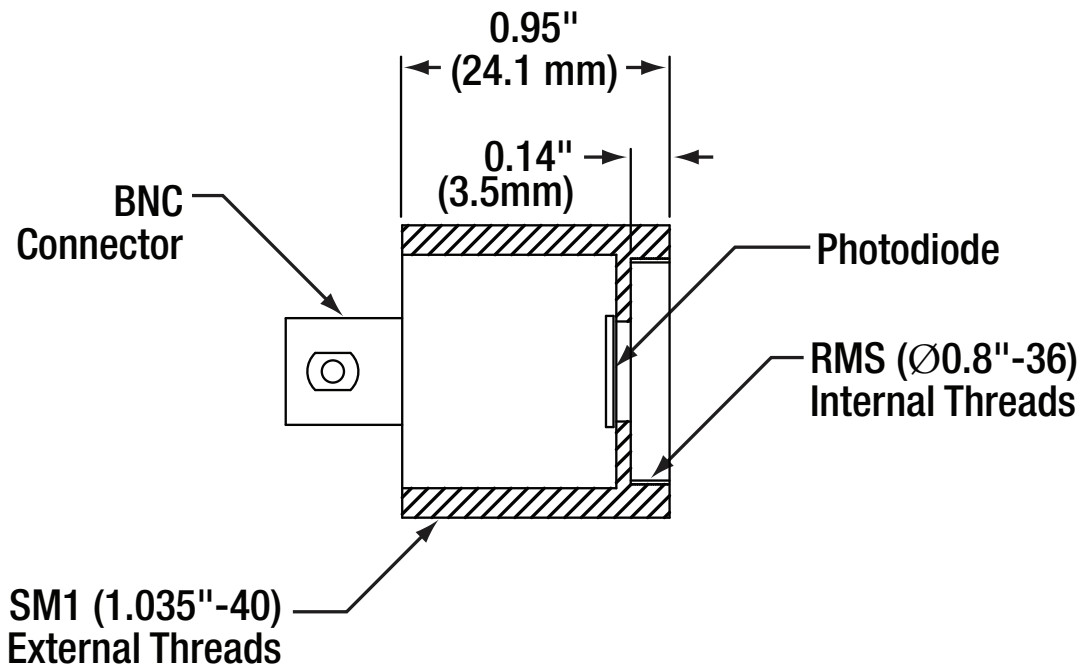

Mounted Photodiode with SM1 Threading

Precautions and Warranty Information

These products are ESD (electro static discharge) sensitive and as a result are not covered under warranty. In order to ensure the proper functioning of a photodiode care must be given to maintain the highest standards of compliance to the maximum electrical specifications when handling such devices. The photodiodes are particularly sensitive to any value that exceeds the absolute maximum ratings of the product. Any applied voltage in excess of the maximum specification will cause damage and possible complete failure to the product. The user must use handling procedures that prevent any electro static discharges or other voltage surges when handling or using these devices.

Thorlabs, Inc. Life Support and Military Use Application Policy is stated below:

THORLABS' PRODUCTS ARE NOT AUTHORIZED FOR USE AS CRITICAL COMPONENTS IN LIFE SUPPORT DEVICES OR SYSTEMS OR IN ANY MILITARY APPLICATION WITHOUT THE EXPRESS WRITTEN APPROVAL OF THE PRESIDENT OF THORLABS, INC. As used herein:

- 1. Life support devices or systems are devices or systems which, (a) are intended for surgical implant into the body, or (b) support or sustain life, and whose failure to perform, when properly used in accordance with instructions for use provided in the labeling, can be reasonably expected to result in a significant injury to the user.*
- 2. A critical component is any component in a life support device or system whose failure to perform can be reasonably expected to cause the failure of the life support device or system or to affect its safety or effectiveness.*
- 3. The Thorlabs products described in this document are not intended nor warranted for usage in Military Applications.*